




WP WP•LS

WessexPower
Solutions for High-tech Instrumentation
www.wessexpower.co.uk

LSTechnology
E-commerce Instrumentation
www.loggershop.co.uk

Telephone +44(0)1929 459 459 Email sales@WPLS.co.uk

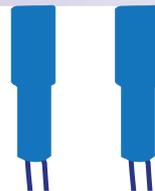

LogTag[®] UTRED30-16

Quick Start Guide

Version H - February 2023


www.logtag.com

Quick Start Guide Contents


What's Included	3
Battery Installation	4
Downloading LogTag Analyzer	6
Configuring your UTRED30-16	7
Starting your UTRED30-16	8
During Recording	10
Downloading Results	11

What's Included


Please check you have each of the items shown below before continuing with setting up your UTRED30-16.

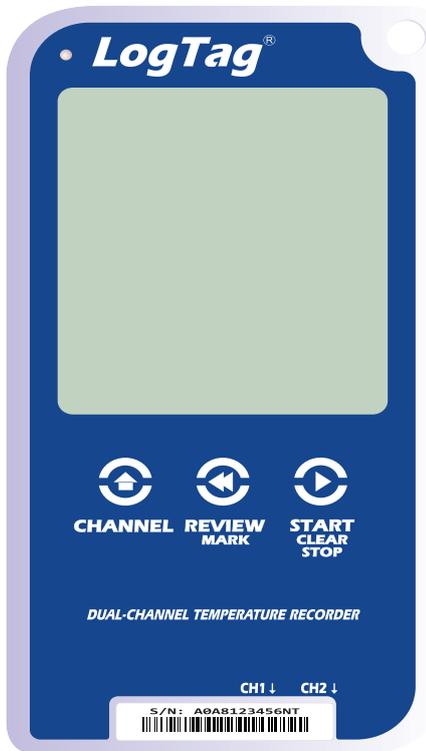

UTRED30-16

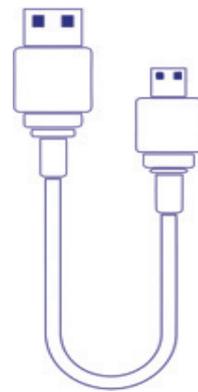


USB Cable

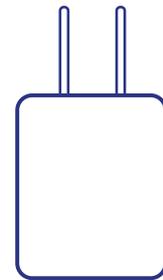


AC Adapter
(US and EU
adapter only)



External Probe
(Not included)

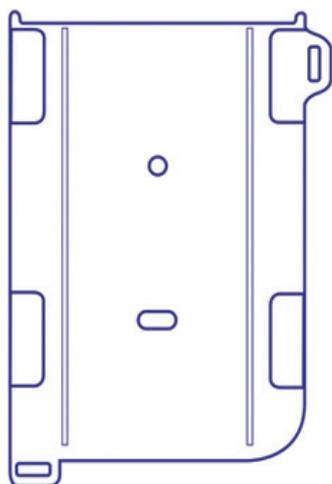

UTRED30
Wall mount
(Not included)

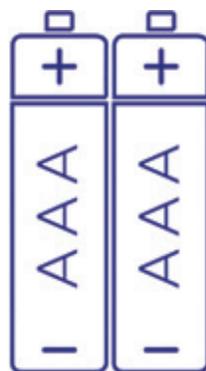

2x AAA Battery
(Not included)



Sensor
Cover
(Optional
to Use)



Glycol Buffer
(Not included)


Battery Installation

The AAA batteries are the main power source for your UTRED30-16.



A USB cable is another power backup source if you wish to power the unit permanently via the USB socket.

1. You will need a small phillips (cross-shaped) screwdriver to remove the cover on the back of the UTRED30-16 case.

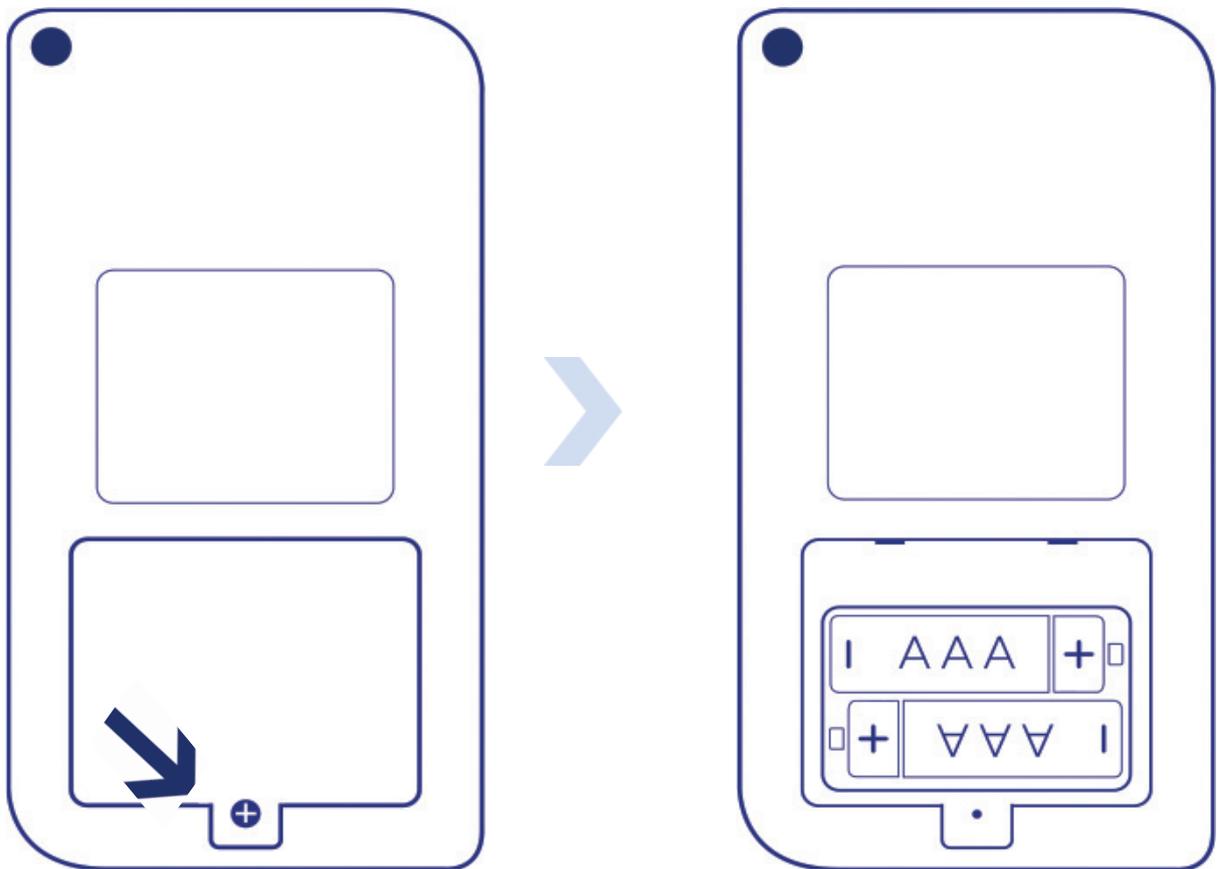

Battery Installation continued...

2. Once you have removed the battery cover, install the 2x AAA batteries.

Take note of the direction each battery must be installed.

3. Once both batteries are securely installed, **replace the battery cover** and secure via the screw you used to open the cover.

Warning: Please note the UTRED30-16 will not start if AAA batteries are low and there is no USB power.

Downloading LogTag Analyzer

To download the latest LogTag Analyzer, open your browser and navigate to:

↓ <https://logtagrecorders.com/software/LTA3/>

1. Click 'Go To Downloads Page' button
2. Click 'Download Now' to start the download
3. Click 'Run' or 'Save File' then double-click the downloaded file to open Analyzer Setup Wizard
4. Once the Analyzer Setup Wizard is open, follow the on-screen prompts to install Analyzer onto your computer
5. Click 'Finish' to exit the Setup Wizard

Warning: Please ensure no other LogTag software is currently running on your computer before running the Analyzer software.

Configuring your UTRED30-16

Connect your UTRED30-16 to your computer via the USB cable provided. The USB socket on the device is located on the bottom, protected by a rubber seal.



1. Open LogTag Analyzer
2. Click 'Configure' from the 'LogTag' menu or click the 'Wizard' icon. Please select which Channel(s) you wish to log, in 'Log channels:' setting.
3. Adjust your logger configuration settings as required. For more information on configuration settings, please refer to [Configuring the UTRED30-16](#) in the [Product User Guide](#) or press 'F1' for help from your keyboard
4. Click 'Next' to upload the configuration settings to the logger
5. Click 'Close' to exit the configuration page

Note: Please ensure you have LogTag Analyzer version 3.1.15 onwards for configuration of the UTRED30-16 dual-sensor models.

Starting your UTRED30-16

Before starting, please ensure you have connected the ST100X external probe(s) into the correct Channel sensor port(s) you have configured to record from.



Press and hold the **START/Clear/Stop** button. **STARTING** will appear along with **READY**.

Release the button once **READY** disappears.

The UTRED30-16 now records temperature data.

The Logger will **not** start if:

- You release the button before **READY** disappears.

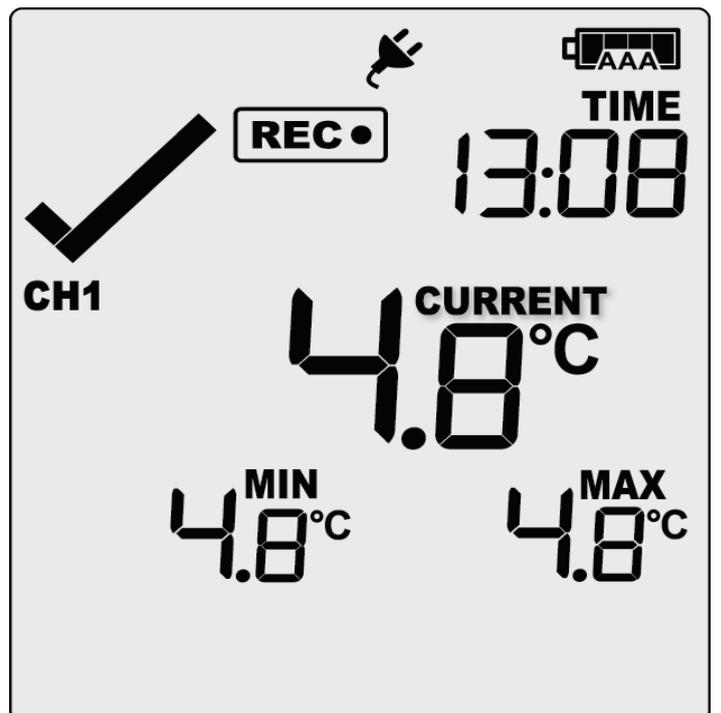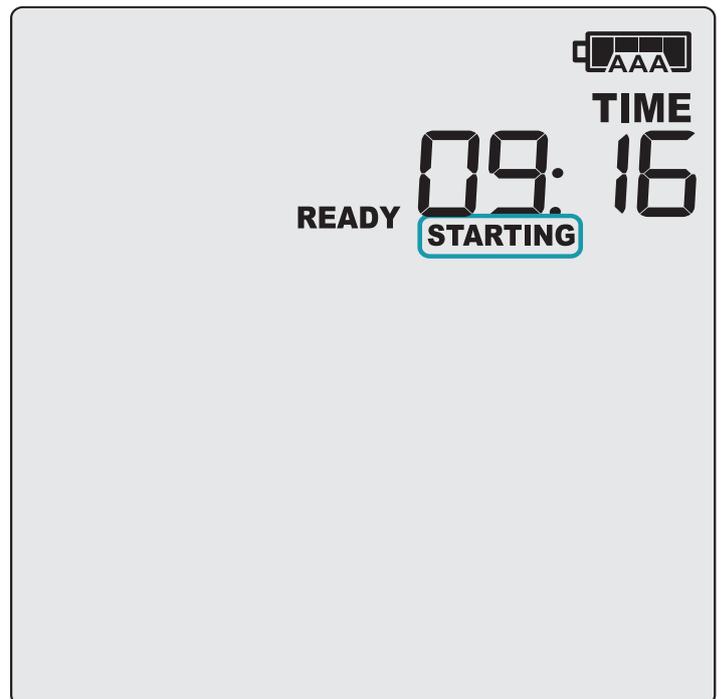

Starting your UTRED30-16 (continued...)

- You keep holding the button for more than 2 seconds after READY disappears.
- The backup battery is critically low and the logger is not connected to power.

Press the **CHANNEL** button to change the channel on screen. In this example, the display switched from CH1 to CH2.



Note: The sensor cover can be inserted in an unused sensor port.

During Recording

To view more information about clearing alarms and resetting and reviewing trip statistics, please refer to the [Product User Guide](#).

End of page

Downloading Results

Connect your UTRED30-16 to your computer via the USB cable provided. The USB socket on the device is located on the bottom, protected by a rubber seal.



A new device drive will appear in file explorer with the files containing the recorded data for the Channel(s) monitored.

Alternatively, you can follow the steps below:

1. Open LogTag Analyzer
2. Click 'LogTag' from the 'Download' menu or press 'F4'
3. Click 'Close' to exit the download page

By default, auto-download is enabled in LogTag Analyzer so when you open the program, data results will be displayed after connecting your UTRED30-16 to your computer